


HPLC DATA SHEET

Sodium Benzoate

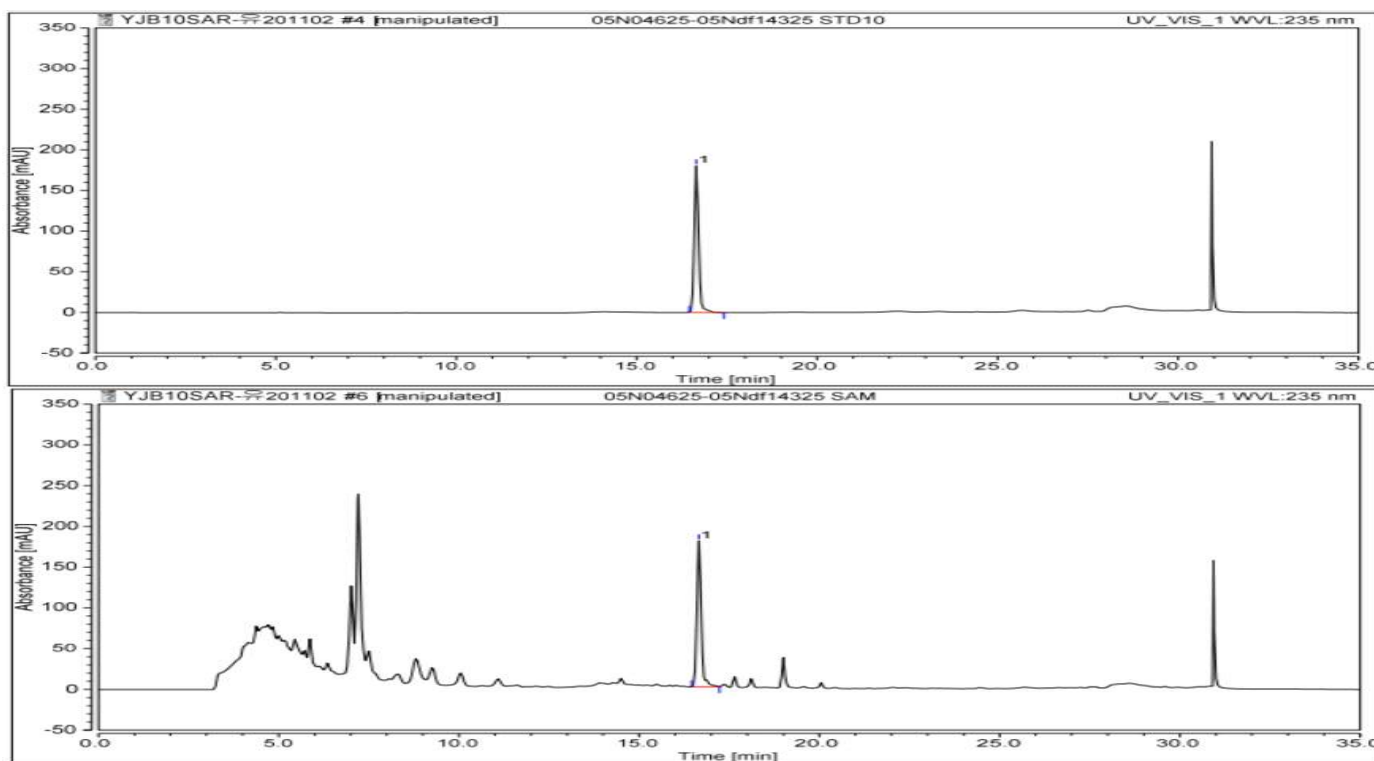

1. Benzoate sodium

HPLC condition

Column INNO Column C18 5um, 4.6mm I.d. × 250mm (P/N 05N04625)

<Gradient system>

	Time (min)	Solution A	Solution B
Mobile Phase	0	100	0
	5	100	0
	25	0	100
	25.1	100	0
	35	100	0

*Solution A : 0.05 % phosphoric acid in 30 % Acetonitrile

*Solution B : 0.05 % phosphoric acid in 90 % Acetonitrile

Flow rate 0.5 mL/min

Temperature 30°C

Detector UV (235 nm)

Injection Volume 20 uL